


“Press-to-Fit” Encapsulating Instructions

Products Shown:

- (1) Part Number: 5415426.1, 5.4mm Sample Pan w/ Flange
- (1) Part Number: 6415426.1, 5.4mm Sample Pan w/o Flange
- (2) Part Number: 5315013.1, “Press-to-Fit” cover
- (1) Part Number: 16T64675, Hand Press tool

Please see the following procedure for press fitting our pan style cover and 5.4mm sample pan.

- 1) Place sample in the center of the pan and then the cover “cup up” with a pair of tweezers
- 2) Turning the tweezers sideways, press the cover down flat, even with top of pan.

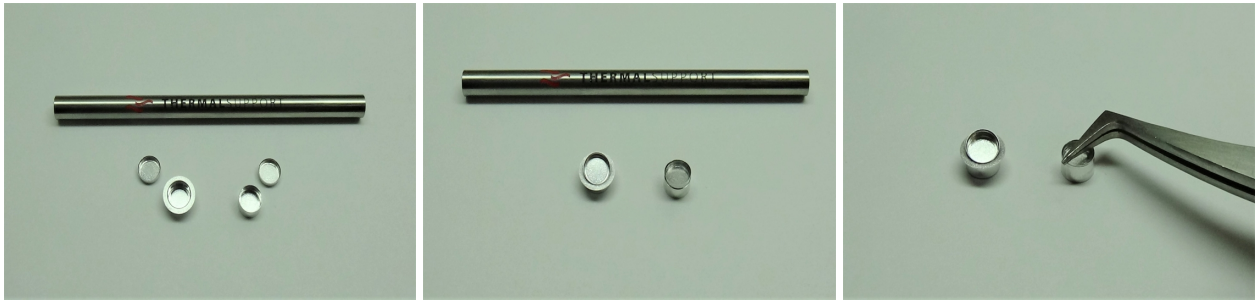

- 3) Using the flat end for films, fibers and powders and the concave end for other samples, firmly press the cover down into the pan with the Hand Press Tool. **Both the pan with flange and the pan without flange are compatible with the TA Tzero crimper with the “Press-to-Fit” cover. Pressing force is stiff but not difficult.*



Please call if you have questions or otherwise need us.