


“Press-to-Fit” Encapsulating Instructions

Products Shown:

- (1) Part Number: 5416717.2, 6.7mm Standard Sample pan
- (1) Part Number: 2416727.2, 6.7mm High Volume Sample pan
- (2) Part Number: 5316500.2, “Press-to-fit” cover
- (1) Part Number: 16T65775, Hand Press tool

Please see the following procedure for press fitting our pan style cover and 6.7mm sample pan.

- 1) Place sample in the center of the pan and then the cover “cup up” with a pair of tweezers
- 2) Turning the tweezers sideways, press the cover down flat, even with top of pan

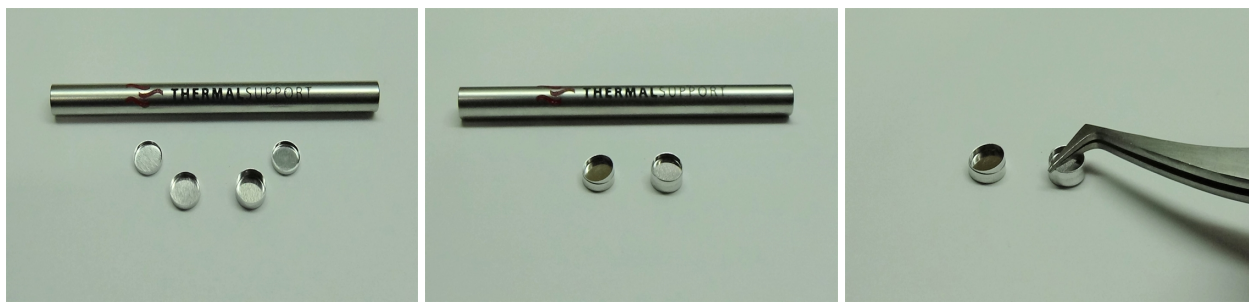

- 3) Using the flat end for films, fibers and powders and the concave end for other samples, gently press the cover down into the pan with the Hand Press Tool. The cover is soft and will typically form around your sample so there is no need to press hard.

Note pushing too hard can damage the bottom of the pan*



Please call if you have questions or otherwise need us.